

MOUNTING INSTRUCTION

Ceiling cover plate for mounting the FX Arm through an inner ceiling

No. 148872/00
Date 2006-12-01

1. Remove a required number of inner ceiling plates.
2. Fasten the bracket against the upper ceiling. Ensure that the surface, on which to mount the bracket is level and use the bracket to mark the holes. Use vibration secured fixing bolts and nuts suitable for the ceiling material. The bolts must each stand a torque force of minimum 7500 N (1650 lbs).
3. If necessary, cut the extension arm to a suitable length.
4. Fit the extension arm to the bracket.
5. Make a hole in the affected inner ceiling plate. Hole diameter: 200 mm. **N.B. The ceiling plate must not be devided.**
6. Refit the lower ceiling plate.
7. Choose a stamped disc on the ceiling cover plate suiting the ventilation duct. Fold it to and fro until it comes loose. Support the discs that are not required to be removed.
8. Fasten the ceiling cover plate. Press it against the inner ceiling.
9. Mount a suitable ventilation duct.
10. Fit a connector to the swivel.
11. Connect the ventilation duct to the swivel.
12. Mount the FX Arm.

